



The easiest protein electrophoresis for increased sample throughput applications



The E-PAGE™ 48 System revolutionizes protein separation and allows rapid, reproducible protein electrophoresis for medium to high sample throughput applications. With E-PAGE™ 48 System, you get:

- **Ease of use**—quick setup and separation of 48 samples or more in just 23 minutes
- **Fast loading**—compatible with multi-channel or robot loading
- **Efficient western blotting and staining**—optimized protocols and reagents

Easy to use, easy to expand

E-PAGE™ 48 8% Gels are self-contained, horizontal pre-cast gels that provide 48 sample lanes and 4 marker lanes with a 3.2 cm separation length. E-PAGE™ 48 Gels are specifically designed for use in E-Base™ devices, which combine an electrophoresis base with a power source. To use, simply insert a gel into a Mother E-Base™ device, load your samples, and run (Figure 1). For higher throughput, you can connect up to three Daughter E-Bases™ to each Mother E-Base™ device (Figure 2) to run a total of 192 samples. Now you can get more results in fewer runs, saving time and effort.

Figure 1 — The easy-to-use E-PAGE™ 48 System


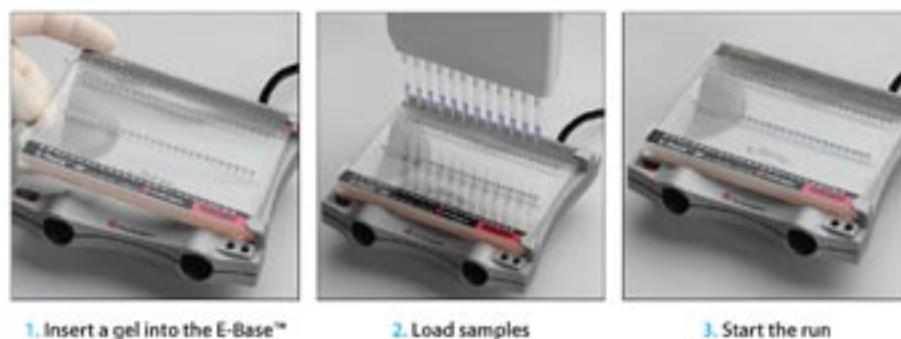


Figure 2 - Easily expand sample throughput using Daughter E-Base™ Devices

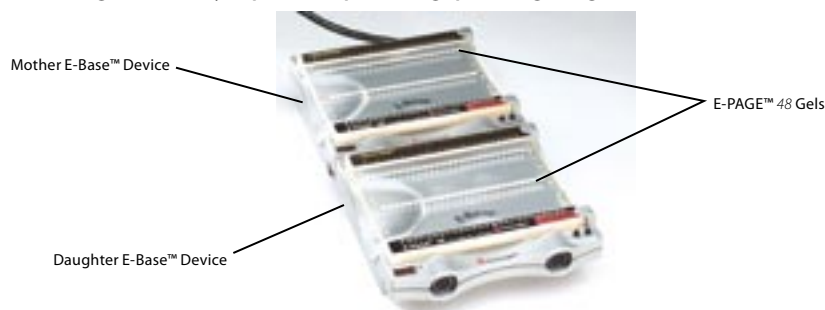

Convenient reagents and software ensure post-run analysis

The E-PAGE™ 48 8% Gel offers a broad molecular weight separation range (Table 1). It's suitable for a variety of post-electrophoresis procedures such as western blotting (Figure 3) and protein staining (Figure 4). Pre-cut blotting membranes, blotting pads, filter papers, blotting roller, and incubation trays (Figure 5) are available to ensure optimal blotting results*. Optimized protocols for Coomassie® and fluorescent staining of proteins are provided to enhance E-PAGE™ staining results. The E-Editor™ v.2 software enables convenient sample lane grouping and alignment for easy band comparison and result interpretation. With the E-PAGE™ 48 System, high-throughput protein analysis has never been easier or more efficient.

Table 1 — Specifications of the E-PAGE™ 48 8% Gel

| | E-PAGE 48 8% Gels |
|----------------------------|--------------------------------------|
| Cassette size | 13.5 (l) x 10.8 (h) x 0.67 cm (w) |
| Gel thickness | 3.7 mm |
| Gel volume | 50 ml |
| Well depth | 3 mm |
| Number of wells | 48 + 4 markers |
| Well format | 2 x 26 |
| Well opening | 3.6 mm x 2.2 mm |
| Running distance | 32 mm |
| Run time | 23 minutes |
| Space between well centers | 4.5 mm |
| Max sample volume | 15 µl |
| Max protein load | 20 µg |

Migration pattern of MagicMark™ XP Standard on an E-PAGE™ 48 8% Gel



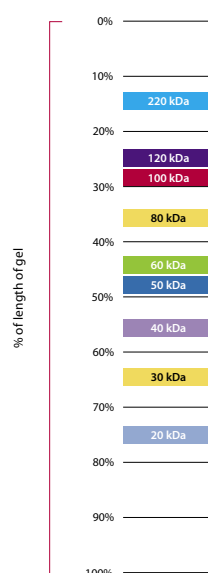


Figure 3 – Western blot results on E-PAGE™ 48 8% Gels

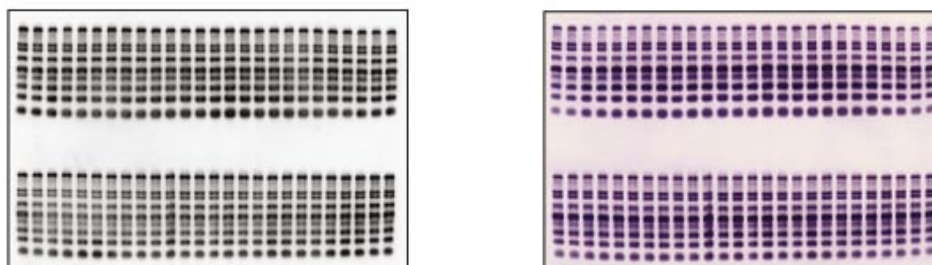

MagicMark™ XP Western Standard (5µl) was loaded into each well of an E-PAGE™ 48 8% Gel and transferred to a pre-cut 0.2 µm nitrocellulose membrane (Cat. no. LC2009). The blot was probed with 1:1000 Anti-Xpress™ Antibody and detected using the WesternBreeze® Immunodetection Kit (anti-mouse) – Chemiluminescent (A) and Chromogenic (B) detection.

Figure 4 — Staining results of E-PAGE™ 48 8% Gel

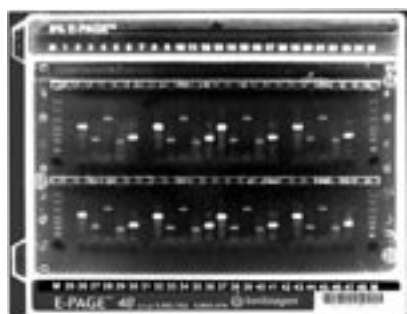

M: BenchMark™ Fluorescent Standard

Lanes 2, 8, 13, 19, 26, 32, 37, 43:

Human kinase fusion protein (10 µl)

Lanes: 3, 9, 14, 20, 27, 33, 38, 44:

E. coli CAT fusion protein (10 µl)

Lanes: 4, 10, 15, 21, 28, 34, 39, 45:

E. coli GUS fusion protein (10 µl)

Lanes: 5, 11, 16, 22, 29, 35, 40, 46:

E. coli calmodulin fusion protein (10 µl)

Lanes: 6, 12, 17, 23, 30, 36, 41, 47:

E. coli kinase D fusion protein (10 µl)

The Lumio-tagged expressed proteins were separated on an E-PAGE™ 48 8% Gel and the entire gel cassette was visualized with a UV transilluminator using an ethidium bromide filter following the recommended protocol in the manual.



Lanes M: 5µl SeeBlue® Plus2 Pre-stained Standard

Lanes 2, 4, 21, 23: 10µl MagicMark™ XP Standard

Lanes 6 & 19: 2µl BenchMark™ His-tagged Standard

Lane 8: Lysozyme 200 ng


Lane 10: Carbonic anhydrase 200 ng

Lane 12: BSA 100 ng

Lanes 15 & 17: 5µl *E. coli* lysate

The gel was stained with Coomassie® R-250 following the recommended protocol in the E-PAGE™ 48 manual.

Figure 5 — E-PAGE™ 48 System and blotting products

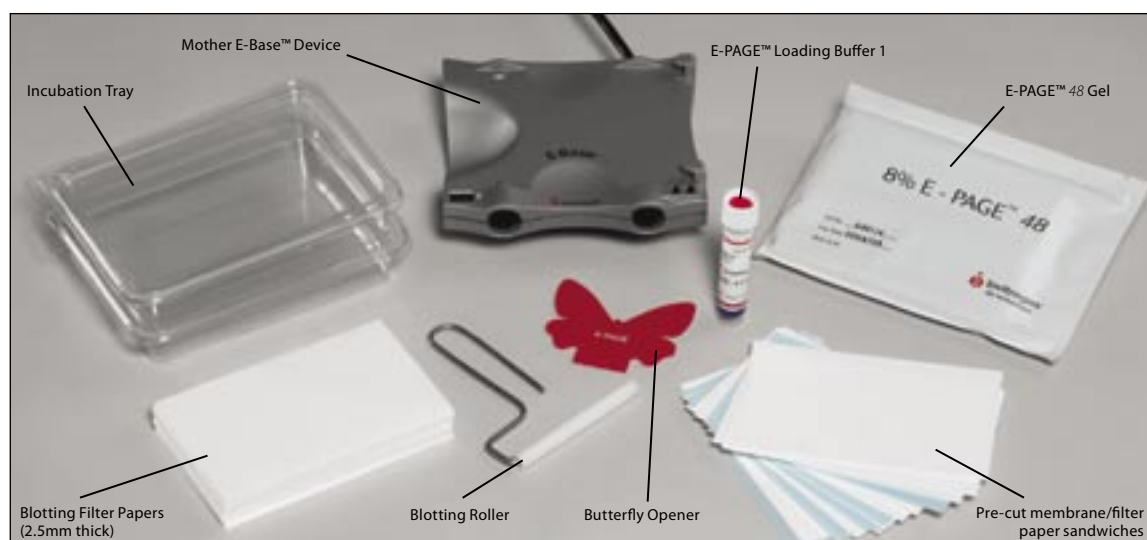

Ordering information

For more information visit www.invitrogen.com/epage. Order the E-PAGE™ 48 System today.

| Product | Quantity | Cat. no. |
|---|---|-----------|
| E-PAGE™ 48 8% Starter Kit | 1 kit | EPST48-08 |
| <i>includes 4 gels, one Mother E-Base™, Loading Buffer 1, one E-PAGE™ Blotting Pad, a Butterfly Opener, and SeeBlue® Plus2 Pre-Stained Standard</i> | | |
| E-PAGE™ 48 8% gels | 8 gels | EP048-08 |
| <i>includes 8 gels, Loading Buffer 1, one E-PAGE™ Blotting Pad, and a Butterfly Opener</i> | | |
| Mother E-Base™ device | 1 unit | EB-M03 |
| Daughter E-Base™ device | 1 unit | EB-D03 |
| E-Holder™ platform | 2 units | EH-03 |
| E-PAGE™ Loading Buffer 1 | 4.5 ml | EPBUF-01 |
| NuPAGE® Sample Reducing Agent (10x) | 250 µl | NP0004 |
| | 10 ml | NP0009 |
| SeeBlue® Plus2 Pre-Stained Standard | 500 µl | LC5925 |
| MagicMark™ XP Western Standard | 250 µl | LC5602 |
| Nitrocellulose/Filter Paper, 8.5 x 13.5 cm size, 0.2 µm pore | 16/pk | LC2009 |
| Nitrocellulose/Filter Paper, 8.5 x 13.5 cm size, 0.45 µm pore | 16/pk | LC2006 |
| Invitrolon™ PVDF/Filter Paper, 8.5x 13.5 cm size, 0.45 µm pore | 16/pk | LC2007 |
| Blotting Filter Paper, 2.5 mm thick | 50/pk | LC2008 |
| Blotting Roller, 8.6 cm wide | 1 unit | LC2100 |
| Incubation Tray, 10 x 14 x 3 cm | 8 trays and lids/pk | LC2102 |
| NuPAGE® Transfer Buffer (20X) | 1 L | NP0006-1 |
| NuPAGE® Antioxidant | 15 ml | NP0005 |
| E-PAGE™ Blotting Pad | 4 pads | LC2101 |
| E-Editor™ 2.0 Software | FREE visit www.invitrogen.com/epage to download | |

Related Products:

WesternBreeze® Immunodetection Chemiluminescent Kits

| | | |
|-------------|-------|--------|
| Anti-Mouse | 1 kit | WB7104 |
| Anti-Rabbit | 1 kit | WB7106 |

WesternBreeze® Immunodetection Chromogenic Kits

| | | |
|-------------|-------|--------|
| Anti-Mouse | 1 kit | WB7103 |
| Anti-Rabbit | 1 kit | WB7105 |

| | | |
|---|-----------|----------|
| Expressway™ HTP Cell-Free <i>E. coli</i> Expression Kit | 5x96 rxns | K9900-80 |
|---|-----------|----------|

| | | |
|----------------------------|-------|--------|
| Lumio™ Green Detection Kit | 1 kit | LC6090 |
|----------------------------|-------|--------|

| | | |
|---------------------------------|--------|--------|
| BenchMark™ Fluorescent Standard | 125 µl | LC5928 |
|---------------------------------|--------|--------|

* Note: a semi-dry transfer apparatus is required to blot an intact E-PAGE™ Gel.

These products may be covered by one or more Limited Use Label Licenses (See the Invitrogen catalog or www.invitrogen.com).
By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses.



For research use only. Not intended for any animal or human therapeutic or diagnostic use.
©2005 Invitrogen Corporation. All rights reserved. Reproduction forbidden without permission. Printed in the U.S.A.

Corporate headquarters:

1600 Faraday Avenue • Carlsbad, CA 92008 USA • Tel: 760 603 7200 • Fax: 760 602 6500 • Toll Free Tel: 800 955 6288 • E-mail: tech_service@invitrogen.com • www.invitrogen.com

European headquarters:

Invitrogen Ltd • Inchinnan Business Park • 3 Fountain Drive • Paisley PA4 9RF, UK • Tel: +44 (0) 141 814 6100 • Fax: +44 (0) 141 814 6260 • E-mail: eurotech@invitrogen.com