

FocalCheck™ Fluorescent Microsphere Standards

Pub. no. MAN0006898

Rev. date 1 October 2012

Detailed protocol is available online at www.lifetechnologies.com/manuals.

WARNING! Before handling, read all applicable Safety Data Sheets (SDSs) at www.lifetechnologies.com/support.

FocalCheck™ Microsphere suspensions

1. Apply 5 μ L of the FocalCheck™ bead suspension to the surface of a slide and spread it with the pipette tip.
2. Wait for the droplet to dry. Apply mounting medium and cover the sample with a coverslip.

FocalCheck™ Microspheres mounted on slides

- One slide contains 6 μ m beads with a green or orange ring stain (Component A).
- Two control slides contain microspheres stained throughout with each green or orange dye separately (Components B and C).

For research use only. Not for use in diagnostic procedures.

Cat. no.	Bead size	Ring stain	General stain
F-7234	15 μ m	blue, orange	—
F-7240	15 μ m	green, dark red	—
F-7235*	15 μ m	green, orange, dark red	—
F-14806**	6 μ m	green, orange, dark red	—
F-7237*	15 μ m	green	blue
F-14808**	6 μ m	green	blue
F-7238	15 μ m	green	dark red
F-7236	15 μ m	orange	blue
F-7239*	15 μ m	dark red	green
F-14807**	6 μ m	dark red	green
F-36905	6 μ m	green-1 (500/512)	green-2 (512/525)
F-36906	6 μ m	orange-1 (532/552)	orange-2 (545/565)

* Also available mounted on slides in Kit F-24634.

** Also available mounted on slides in Kit F-24633.

† Only available mounted on slides as a kit.

Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

© 2012 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

DISCLAIMER — LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

For support visit www.lifetechnologies.com/support

www.lifetechnologies.com