

# MultiShot™ StripWell TOP10 Chemically Competent *E. coli*

Catalog Numbers C4096-01, C4096-05, C4096-10

Document Part Number 250374

Publication Number MAN0000192

Revision 2.0

## Description

The MultiShot™ StripWell TOP10 Chemically Competent *E. coli* Kit is chemically competent TOP10 *E. coli* packaged in a 96-stripwell plate to simplify medium- and high- throughput bacterial transformation. The stripwell format allows you to use only the number of wells you need for your particular application.

## Contents

Fifty microliters of chemically competent TOP10 *E. coli* are supplied per well with a transformation efficiency of  $> 1 \times 10^8$  cfu/ $\mu$ g pUC19 DNA. Each MultiShot™ StripWell Kit contains the following reagents.

Catalog no.	Number of plates	S.O.C.	pUC19
C4096-01	1	2 × 15 mL	50 $\mu$ L (500 pg)
C4096-05	5	10 × 15 mL	50 $\mu$ L (500 pg)
C4096-10	10	20 × 15 mL	2 × 50 $\mu$ L (1 ng)

## Genotype

*F- mcrA Δ(mrr-hsdRMS-mcrBC) φ80lacZΔM15 ΔlacX74 recA1 araD139 galU galK Δ(ara-leu)7697 rpsL (Str<sup>R</sup>) endA1 nupG*

## Shipping and Storage

Each MultiShot™ StripWell Kit is shipped on dry ice. Upon receipt, store the kit at  $-85^{\circ}\text{C}$  to  $-68^{\circ}\text{C}$ .

**Note:** You may store the S.O.C. medium at  $15^{\circ}\text{C}$  to  $30^{\circ}\text{C}$  or at  $2^{\circ}\text{C}$  to  $8^{\circ}\text{C}$ .

## Composition of SOC Medium and pUC19

### S.O.C. Medium:

2% Tryptone; 0.5% Yeast Extract; 10 mM NaCl; 2.5 mM KCl; 10 mM MgCl<sub>2</sub>; 10 mM MgSO<sub>4</sub>; 20 mM glucose.

### pUC19:

Each tube contains 10 pg/ $\mu$ L pUC19 in 5 mM Tris-HCl, 0.5 mM EDTA, pH 8.

## IMPORTANT!

Do not transform chemically competent cells by electroporation. The salt content of the buffer will cause arcing and kill the cells.

## Procedure introduction

The following procedure describes how we qualify the MultiShot™ StripWell Kit. In addition, this information will help you determine how to best use the MultiShot™ StripWell Kit for your own applications.

## Total well volume

Each well holds 1.2 mL.

## Before you begin

- Prepare a container of ice large enough to chill the number of wells you will be using.
- Bring the vial of S.O.C. to room temperature.
- Pre-heat a water bath to  $42^{\circ}\text{C}$ .

## Procedure

1. Remove a MultiShot™ StripWell plate from the freezer and remove the number of wells you need. Return any unused wells to the freezer. Place the wells in the container of ice. The cells should thaw within 1 minute.
2. Carefully remove the strip of caps from each set of 8 wells and keep them for further use.
3. Use a multi-channel pipette to add 2–5  $\mu$ L DNA (2 pg–20 ng) to the wells. Keep the volume around 2  $\mu$ L for uniform results.
4. After adding the DNA, cover the wells with the caps and incubate the cells and DNA on ice for 30 minutes.
5. Transfer the wells to the water bath and heat-shock them for 30 seconds at  $42^{\circ}\text{C}$ .  
**Note:** Be careful not to contaminate the cells.
6. Transfer the wells back to the ice and allow the wells to cool for 1 minute.
7. Remove the caps and add 250  $\mu$ L S.O.C. to each well. Re-cap the wells tightly.
8. Incubate the wells at  $37^{\circ}\text{C}$  for 1 hour with shaking (225 rpm). We turn the wells on their side to increase aeration and secure them to the shaker.
9. Plate the appropriate volume from each well. See the next page for examples and expected number of colonies.

## Plating volumes and expected results

The following table describes the type of DNA, amount transformed into MultiShot™ StripWell chemically competent cells, the volume plated, and the number of colonies.

**Note:** We use pUC19 to qualify the kit. Transformation efficiency should be  $> 1 \times 10^8$  cfu/ $\mu\text{g}$  and yield 100–300 colonies per plate. Variability should be no more than 5-fold between wells.

DNA	Type	Amount Transformed	Volume Plated	Number of Colonies
pUC19	Supercoiled	5 pg plasmid	10 $\mu\text{L}$	100–300
pCR™2.1-TOPO® plus amplified 750 bp insert	TOPO® Cloning Vector	3.4 ng vector + insert	25 $\mu\text{L}$	100–300

## Safety Data Sheets (SDSs)

Safety Data Sheets (SDSs) are available at [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support).

## Certificate of Analysis

The Certificate of Analysis provides detailed quality control and product qualification information for each product. Certificates of Analysis are available on our website. Go to [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support) and search for the Certificate of Analysis by product lot number, which is printed on the box.



## Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at [www.lifetechnologies.com/termsandconditions](http://www.lifetechnologies.com/termsandconditions). If you have any questions, please contact Life Technologies at [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support).

## DISCLAIMER

LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.